

B6-8286 ANDROID TABLET CAMERA INSTALLATION INSTRUCTIONS

Procedure:

C-Mount Adapter Location: The microscope is equipped with a pre-mounted C-mount adapter located on the top of your microscope. No adjustments to this adapter are required.

Protective Cap Removal: Remove the protective cap covering the C-mount adapter's threaded opening.



Camera Insertion: Align the camera's threaded connector with the C-mount adapter's threads.

Camera Tightening: Rotate the camera clockwise to engage the threads. Tighten until a snug fit is achieved. Avoid over-tightening. This completes the assembly of the B6-8286 Android Camera Tablet.



UNICO

182 Ridge Road, Suite E, Dayton, NJ
08810 USA Email: info@unicosci.com
Phone: 732-274-1155

unicosci.com