



## Accessory Installation Guide Water-Jacketed Single Cuvette (S2100-105P)

- Step 1: Remove all items from the accessory package made sure that they are complete:
- *A Universal Platform* 1 pc.
  - *Water-Jacketed Holder* 1 pc.
  - *Flow-thru Panel* 1 pc.
  - *Installation Guide* 1 pc.
- Step 2: The only tool needed is a standard “Philips” head screwdriver.
- Step 3: Remove the 4-cell sample holder: Open the sample compartment cover and remove any cuvettes or any breakable items. Push-in the cell holder rod (black knob) all the way back (First position cell holder is into the light path).
- Step 4: Unscrew the cell holder rod by turning the knob counterclockwise until removed from holder. Gently slide the guide rod out of the sample compartment.
- Step 5: Looking over the sample compartment, remove the front screw (Figure 1).

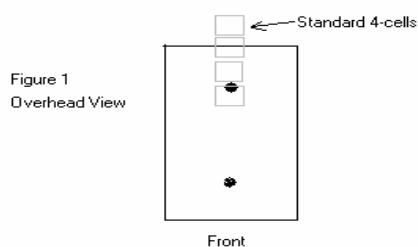


- Step 6: To remove the back screw, place your hand ON TOP of the 4-cells and gently slide towards the front. The back screw is then exposed and can be removed.
- Step 7: Insert Universal Holder by repeating Steps 4-6 in reverses.
- Step 8: Screw in the Water-Jacketed Holder to the **first** (front) position of the Universal Holder (platform). This is to prevent stretching the tubing.
- Step 9: Remove the front panel by sliding the panel up and out (Figure 2).

Figure 2: Removable Front Panel  
Front View

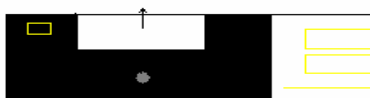

- Step 10: Insert Front Panel for Water-Jacketed Cuvette in place of removed panel.
- Step 11: Connect Water bath to the front panel tubing connectors (Figure 3).

Figure 3: Front Panel for  
Water-Jacketed Cuvette Accessory

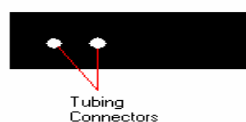

- Step 12: Set desired temperature on water bath (see instructions of water bath manufacturer).