



Eyepiece Digital Camera
Operation Manual
For use with MFR#B6-8182



United Products & Instruments, Inc.
182 Ridge Road Suite E, Dayton NJ USA
(732)274-1155
info@unicosci.com
www.unicosci.com

UNICO Digital Eyepiece Camera Specifications

Sensor Size	1/3.2"
Resolution	1600x1200 (5FPS) 1280x1024 (7.5FPS) 1280x960 (7.5FPS) 800x600 (20FPS)
Exposure	Auto
White Balance	Auto
Power Input	DC 5V (via USB 2.0 PC port)
Eyepiece Size	23mm
Operating System	MS Windows XP/Vista/7/8/10 32 & 64bit

Camera Package includes

Digital Eyepiece Camera	1 piece (body color may vary)
USB Cord	1 piece
Software CD	1 piece

Camera Installation

1. Remove the eyepiece from your microscope and store it in a clean dry place
2. Insert the camera into the eye port of your microscope
3. Connect the USB cord to the camera and your PC
4. Run the software installation from the CD included
5. Start the software and follow user manual included on CD

When finished – remove the camera and replace the eyepiece.

Make sure camera is stored in a clean dry place when not used.

